


# BYO Filter Kit Sampling Instructions




If you would like more information including notes on sampling, please scan the code to visit [wilderlab.co/directions](http://wilderlab.co/directions)



1 Wearing the provided gloves, use the sterile tweezers to place filter into the Falcon tube.

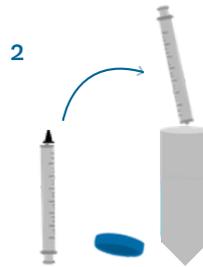

2 Remove the black cap from the small syringe and dispense the entire preservative into the Falcon tube.

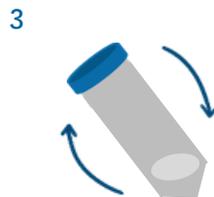

3 Recap the Falcon tube and gently shake to evenly coat the filter with the preservative.



4 Place sealed Falcon tube back into the same kit bag. Ensure the UID label remains visible. **NOTE: Please ensure only 1 filter and associated Falcon tube is placed in each kit bag.**



5 Seal the sample bag and record the sample details in the space provided. Ensure that the coordinates are entered in WGS84 decimal format (for example -41.30951, 174.82110 as displayed on Google Maps).



**Submit your samples online at [wilderlab.co/submit-samples](http://wilderlab.co/submit-samples)**

Print and sign the Chain of Custody (CoC) form that will be emailed to you after online sample submission. Include this in the parcel containing your samples (no refrigeration necessary).

**Send your samples back to us within one month from collection**

If this isn't possible, ensure that your samples are stored in a freezer at -20 C until you send them.

**Send your samples and CoC form to the address provided in the auto-email**