

# eDNA Swab Kit Instructions



If you would like more information including notes on sampling, please scan the code to visit [wilderlab.co/directions](https://wilderlab.co/directions)

1




Take the gloves out of the sample bag and put them on. Pull the swab stick out of the tube and swab the sample.

2



Unscrew the lid of the micro tube and place the swab stick in the preservative.

3

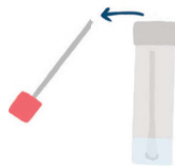

Snap off the end of the swab stick and discard this part. Leave the rest of the swab in the tube and screw on the lid.

4



Place the micro tube into the sample bag. Seal the sample bag and record the sample details in the space provided



**Submit your samples online at [wilderlab.co/submit-samples](https://wilderlab.co/submit-samples)**

Print and sign the Chain of Custody (CoC) form that will be emailed to you after online sample submission. Include this in the parcel containing your samples (no refrigeration necessary).

**Send your samples back to us within one month from collection**

If this isn't possible, ensure that your samples are stored in a freezer at -20 C until you send them.

**Send your samples and CoC form to the address provided in the auto-email**